


BROEN-LAB

recessed eye wash & body shower - mounting instructions

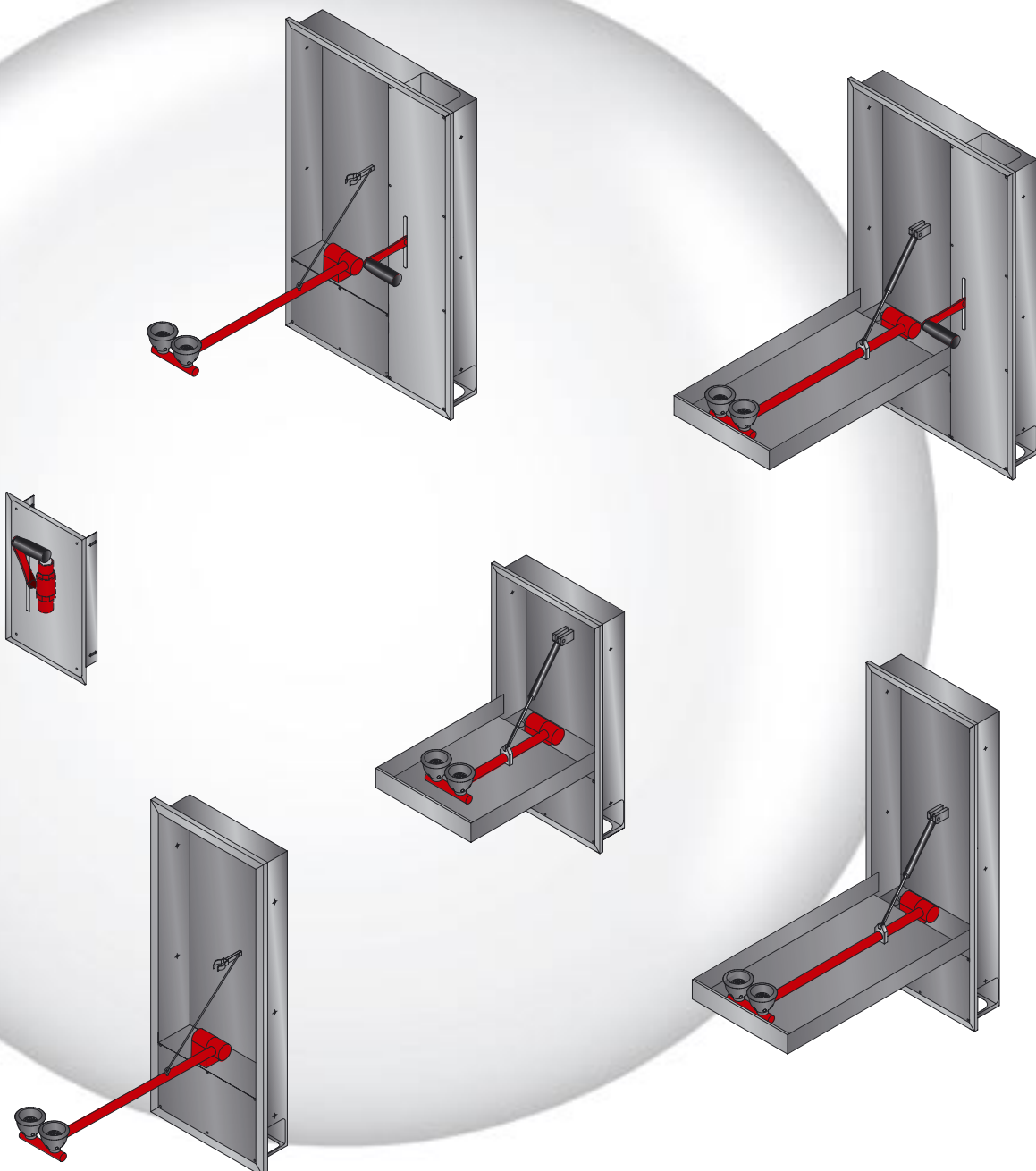


BROEN-LAB

eye washes


- mounting instructions

For models: 17 151.009 / 17 152.009 / 17 161.009 / 17 162.009


Hole in wall, dimensions

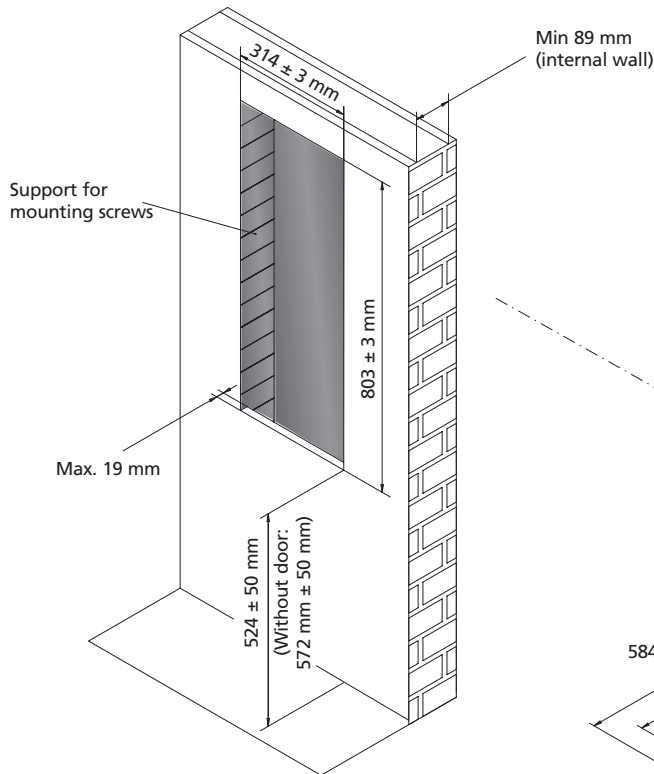

Cabinet, dimensions

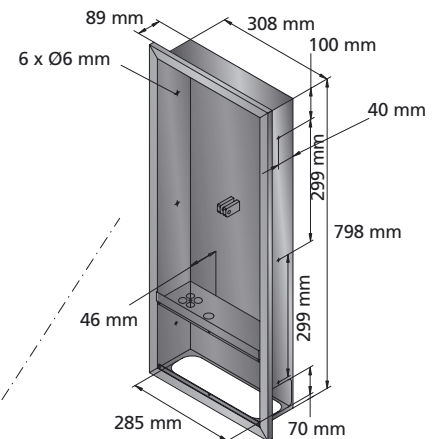


Supply, dimensions

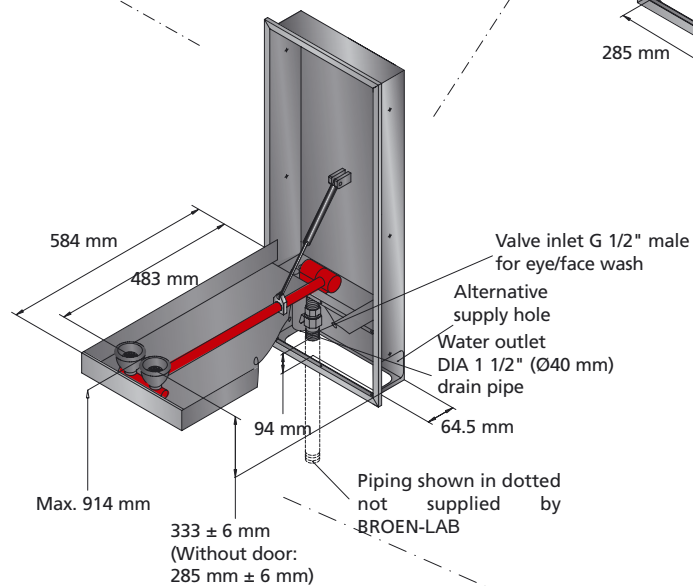


Surface mounting

The cabinet can be surface mounted on the wall by drilling 4 holes in the back wall of the cabinet. Please be aware that the piping will be exposed by surface mounting.

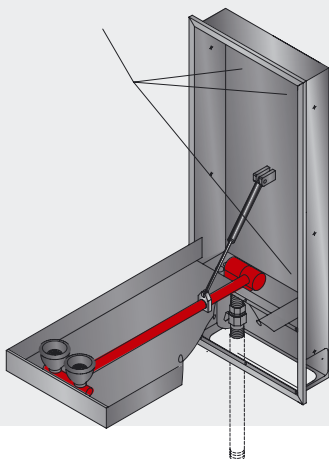

STABITHERM supply, dimensions

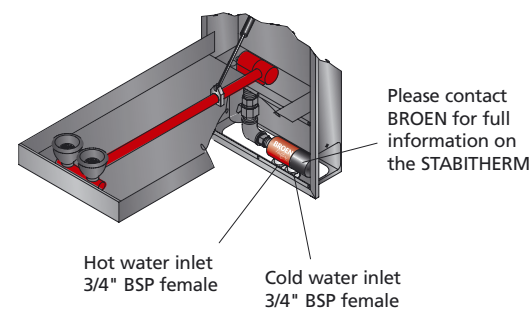


Please contact BROEN for full information on the STABITHERM

- body showers

- mounting instructions

For models: 17 155.009 / 17 165.009

Hole in wall, dimensions

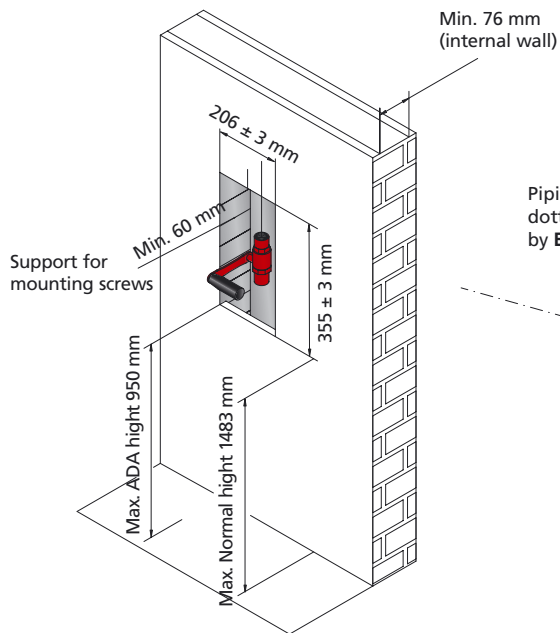

Supply, dimensions

Piping shown in dotted not supplied by **BROEN- LAB**


Valve inlet and outlet G 1" female for body shower

269 ± 5 mm


For ADA purpose Max: 1219 mm

For normal purpose Max: 1752 mm

Cabinet, dimensions

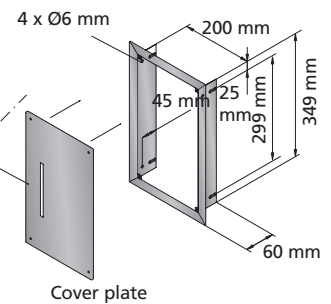


Handle to be mounted after the cover plate

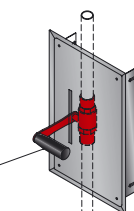

Tepid water outlet G 1" female for body shower






Tepid water part No. 17 457.009

Hot and cold water inlet Pipefix f. Ø28 mm

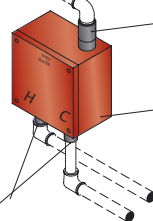

For models: 17 153.009 / 17 154.009 / 17 163.009 / 17 164.009

Hole in wall, dimensions

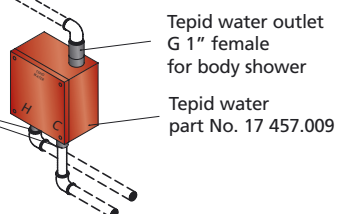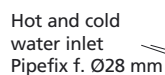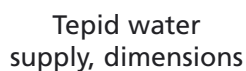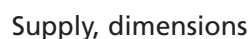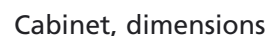



BROEN-LAB manufacture and market a wide range of quality products for high performance flow regulation. The products primarily apply to the professional user in all types of laboratories, the educational sector, the health sector, the pharmaceutical industry and industry in general.



**BROEN-LAB
ISO 9001 CERTIFICATION**
In September 1991 BROEN LAB GROUP was certified according to ISO 9001 as one of the first Danish companies. The certification was carried out by Bureau Veritas, Denmark.

The quality management system of BROEN-LAB now complies with detailed specifications laid down by the internationally acknowledged EN ISO9001:2000. This certification will further contribute to BROEN-LAB's reputation for quality and reliability.

BROEN-LAB

broen-lab.com

HEADQUARTER

BROEN-LAB A/S

Drejervaenget 2 · DK-5610 Assens
Tel. +45 6376 6376 · Fax +45 6471 2476 · broen-lab.com · lab@broen-lab.com

INTERNATIONAL OFFICES

BROEN-LAB Ltd.

Unit 9 Navigation Point · Golds Hill Way · Tripton · West Midlands DY4 OPY · UK Tel.
+44 121 522 4515 · Fax +44 121 522 4535 · broen-lab.com · lab@broen-lab.com

BROEN-LAB Ltd.

10 Bukit Batok Crescent # 10-02 · The Spire · Singapore 658079
Tel. +65 6298 0662 · Fax +65 6298 0468 · broen-lab.com · lab@broen-lab.com

BROEN-LAB GmbH

Tel. +45 6376 6376 · Fax +45 6471 2476 · broen-lab.com · lab@broen-lab.com

BROEN-LAB Inc.

15 Constitution Drive · Suite 122 · NH 03110 · Bedford · USA
Tel. +1 603 310 5089 · broen-lab.com · lab@broen-lab.com

